



Eden Materials

## **Application Note: Flexdym™ Chip Bonding**

Assembly & Bonding of Flexdym™ Microfluidic Devices

Issue Date: 18/04/2020



### **Eden Tech**

Registered office: 83 Avenue Philippe Auguste, Paris 75011, France

Registration / SIRET: 83325407100016

# 1 Summary

This application note describes the bonding of Flexdym™, a proprietary styrenic polymer, developed by Eden Tech. Flexdym™ has mechanical and surface properties which allow it to bond to itself and a variety of other materials, including polystyrene (PS), cyclic olefin copolymer (COC), polycarbonate (PC), poly(methyl methacrylate) (PMMA) and glass. Consequently, Flexdym™ can be useful for various applications ranging from cosmetics and environmental analysis, to organ-on-chip, diagnosis, drug discovery, medical device, etc.

One of the key advantages of using Flexdym™ is its easy and user-friendly bonding method. Unlike most microfabrication materials, it does not involve a time-consuming protocol, expensive material, nor any harsh chemical treatments. There is no oxygen plasma treatment required, in contrast to PDMS, one of the most popular materials currently used for microfluidics prototyping.

The Flexdym™ material bonds to its substrate directly on contact. This procedure may be performed at room temperature, or with heat added to accelerate the bonding process. If it no longer has any use, the Flexdym™ may also be un-bond from its substrate. This recyclable characteristic of Flexdym™ also renders it eco-friendly and low cost. Used Flexdym™ may be sterilized and re-molded to create new microfluidics chips with different designs.

# 2 Introduction

Easy bonding using Flexdym™ is possible thanks to its mechanical and surface properties which have been characterized. The Young modulus of Flexdym™ is 1.15 MPa, encouraging intimate contact with the surface of the substrate it is bond to.

Hardness Shore A	35
Specific Gravity	0.9 g/ cm <sup>3</sup>
Tear Strength	15 kN/ m
Tensile Strength	7.6 MPa
Young Modulus	1.15 Mpa
Elongation	720%
Melt Flow Rate	2 g/ 10 min.

Flexdym™ is a moderately hydrophobic material with dynamic contact angles (CA) measured at:

Advancing CA	105° ± 4°
Receding CA	88° ± 4°

Following oxygen plasma treatment, the static contact angles of Flexdym™ demonstrate stability for up to 4 days (time period limited by experiment length). This shows Flexdym™ has higher hydrophilic stability than PDMS.

Flexdym™ can bind to most substrates at room temperature thanks to its adhesive properties. The bonding strength of Flexdym™ depends on 3 key parameters: time, temperature and surface contact.

Flexdym™ can be strongly bond to itself, and a range of other polymers, including polystyrene (PS) and cyclic olefin copolymer (COC), withstanding fluidic pressures up to 2 Bar. Additionally, it can also be bond to glass substrate with moderate bonding strength.

Interestingly, Flexdym™ may be weakly bond to skin, potentially serving a unique function as a microfluidic skin patch.

For users interested in recycling their Flexdym™ devices, it is important to know that Flexdym™ binds to itself permanently.

Flexdym™ bonding to other polymers is reversible, however it becomes increasingly difficult as the time of contact increases. Reversible binding is facilitated the most by glass substrates.

However, it is important to note this bonding behavior may vary depending on the material grade. If required, the bonding strength can be enhanced by using a coating agent on the substrate.

The Flexdym™ has 2-years shelf life warranty.

### 3 Overview of Flexdym™ Bonding & Assembly

The bonding of Flexdym™ is a straightforward procedure, requiring one step: gently stack the Flexdym™ sheet on the substrate material by manual or automated lamination.



For glass substrates, bonding strength may be enhanced via oxygen plasma.

**Optional :** To accelerate bonding, heat the device using either a hot plate or oven. Then cool down at room temperature.

## 4 Temperature & Time Bonding Parameters

A key advantage of Flexdym™ is its ability to bind substrates at room temperature. This is ideal for biological applications, where damage of reagents or surface modifications must be avoided. Bear in mind, in this case the time required to bond Flexdym™ must be increased.

However, the bonding process can also be accelerated via heat treatment. This simply involves heating the device in an oven or a hotplate. For improved surface contact between Flexdym™ and the substrate, a weight may be required.

The table below shows the bonding parameters required for accelerating Flexdym™ bonding to different substrates:

Substrate	Temperature (°C)	Time (min.)	Bonding Strength
Flexdym	80-85	30	Excellent
Thermoplastics, PS & COC	95	60	Excellent
Thermoplastics PC & PMMA	95	60	Moderate
Glass	95	60-120	Moderate

## 5 References

Lachaux, J., *et al.* (2017) Thermoplastic elastomer with advanced hydrophilization and bonding performances for rapid (30 s) and easy molding of microfluidic devices. *Lab on a Chip* 17, 2581-2594.

Roy E., *et al.* (2015) From cellular lysis to microarray detection, an integrated thermoplastic elastomer (TPE) point of care Lab on a Disc. *Lab on a Chip* 15, 406



Eden Materials is a division of Eden Tech  
Registered Address: 83 Avenue Philippe Auguste, 75011 Paris, France  
Client Facing Address: 172 Rue de Charonne, 75011 Paris, France  
T; + (33) 188 327 405  
E; [contact@eden-microfluidics.com](mailto:contact@eden-microfluidics.com)  
W: [www.eden-microfluidics.com](http://www.eden-microfluidics.com)